

How to Change the Tamp Pad on a Sci-Print MP2/MP2+

Purpose


This guide shows the proper way to change the tamp pad for the Sci-Print MP2/MP2+. To purchase the Standard Tamp Pad Kit (1043-310-101) that is used to replace a damaged tamp pad, please contact sales@scinomix.com or 314-298-9800 Option 1. To replace the damaged tamp pad, please follow the steps listed below. Should you need further assistance, please contact service at 314-298-9800 ext. 208 or e-mail service@scinomix.com.

Who Should Use This Troubleshooting Guide?

End User

Instructions

1. Turn off air supply to the unit.
2. Locate supplied replacement Tamp Pad Kit (part number 1043-310-101). The Standard Tamp Pad Kit includes 2 Standard Tamp Pads, 0.25" x 2" (part number 1043-310-011), 2 Tamp Pad Covers (part number 1043-310-013), and 3 Positioning Pins (part number 2000-128).

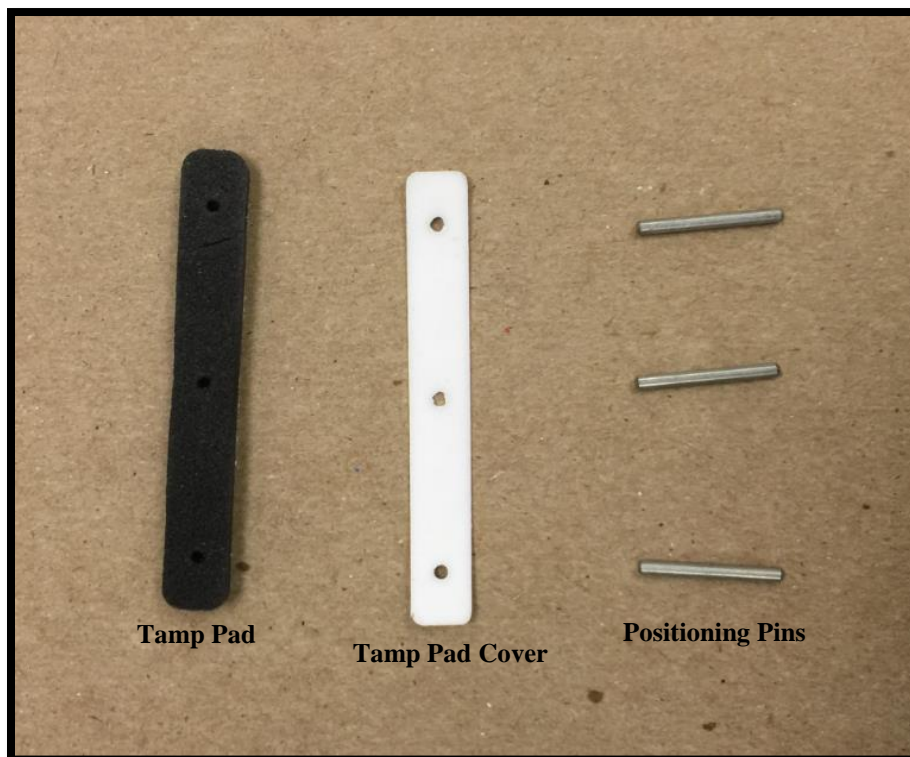

3. Remove the Tamp Pad from the tamp head using fingers or a flat head screw driver. (Be careful not to damage the tamp head, and only remove the foam).



4. Wipe off any excess glue with an alcohol pad leaving the tamp head completely clean.



5. Put positioning pins provided through the 3 holes on the Tamp Pad, but do not remove the white backing covering the adhesive.




6. Remove backing from Tamp Pad Cover.
7. Place Tamp Pad Cover over the Positioning Pins, with the adhesive side facing the non-adhesive side of the foam, and seat onto the Tamp Pad.



8. Remove backing of Tamp Pad.

9. Place Positioning Pins into holes of tamp head

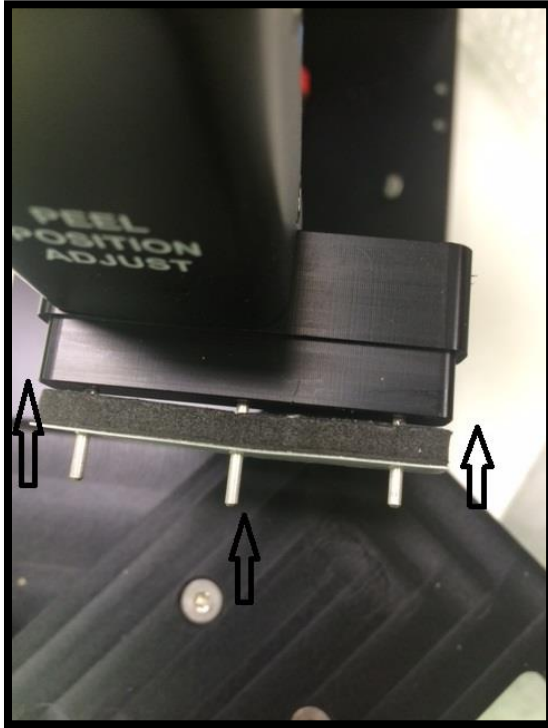

10. Push the pad onto the tamp head.
11. Remove the Positioning Pins



12. Turn air to unit on, and test using the diagnostic screen.

TS-1027
10/2015



Version Number	Change	Date Effective	Approved by:
A	New	10/2015	EP